

# File transfer from analysis cluster

Editors: Changwoo Do, Gergely Nagy, Yingrui Shang, William Heller  
Date: 07/26/2021

## Introduction

After reducing the data files with any of the following methods

- via scripts run in the terminal window,
- via mantidworkbench or
- via jupyter notebooks

the reduced files can be transferred from the analysis cluster to your personal computer. The content of the document is also explained through a tutorial video: <https://www.youtube.com/watch?v=e01ac0Nz6DQ>

Below please find the step-by-step description of the procedure.

### 1. Visit [analysis.sns.gov](https://analysis.sns.gov)

- Log in with your xcams ID and password (IPTS system username and password)


### 2. Zip and send the selected data files

- Open a terminal
- Access the folder where the reduce files were saved with the use of:  

```
cd /SNS/EQSANS/IPTS-yourIPTNumber/shared
```
- Data can be zipped with the zip command. E.g. zipping all the files containing the word merged can be done via:
  - zip data.zip merged\*
- The zipped file can be mailed to yourself with the title "data from IPTS-**yourIPTNumber**" via:
  - echo "data from IPTS-**yourIPTNumber**" | mail -s "title" -a data.zip youremail@email
  - mail -s "title" -a data.zip youremail@email <<< 'This is the email text'

### 3. Alternative solution. Use the freely available WinSCP client

- Download the WinSCP client from <https://winscp.net/eng/index.php>
- Install the software
- Open the program, you will see a login window as below

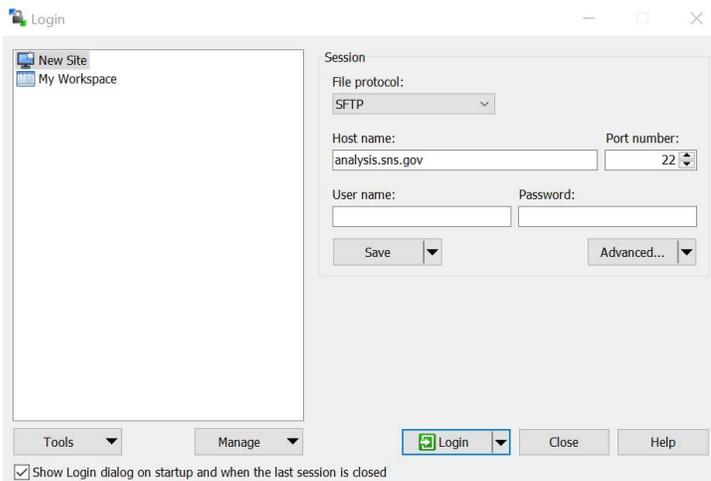

- For hostname use `analysis.sns.gov`
- For username and password use your xcams ID and password
- Access `/SNS/snfs1/instruments/EQSANS/IPTS- yourIPTNumber/shared`

- g. Access the folder selected during data reduction and copy the files of interest to your computer
- h. The + button above the analysis cluster side of the window will allow you to select files containing a specific textSegment via typing \*textSegment\* into the window

**4. FileZilla is another popular software for FTP/SFTP file transfers**

- a. It can be downloaded from <https://filezilla-project.org/>
- b. To connect to analysis cluster
  - i. Click on Edit->Site manager, on the left panel click on new site
  - ii. On the right panel:
    1. Protocol: SFTP
    2. Host: analysis.sns.gov in Host
    3. User & Password: xcams username and password
    4. Connect