
HIO-QSG-023 Page 1 of 1

Super-Nuova Milti-Place Stirrer and Stirring Hot Plate

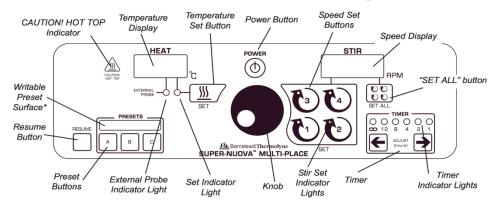



Temperature range: 1°C − 370°C; Stirring speed range 50 − 1200rpm;

General Operation

- 1. Turn on the power button which is located on the front control panel;
- 2. **To set the individual stirring speed**, press desired stirring position key (1-4) located under the "STIR" display. Use the knob in the center of the console to select a speed, turn the knob clockwise to increase the speed or counterclockwise to decrease the speed. When desired speed has been reached, press same key (1-4) under the "STIR" display or wait a moment and the unit will beep indicating that it has been set. Repeat for each individual stirring position.
- 3. **To set all positions to the same stirring speed**, press the "SET ALL" key under the "STIR" display, use the knob in the center of the console to select a speed. Turn the knob clockwise to increase the speed or counterclockwise to decrease the speed. When desired speed has been reached, press the "SET ALL" key again or wait a moment and the unit will beep indicating that it has been set.
- 4. **To set the temperature**, press "SET" key located under the Heat display. The red set temperature indicator light will blink, indicating it is ready to be set, use the knob in the center of the console to select a temperature. Turn the knob clockwise to increase the temperature or counterclockwise to decrease the temperature. When desired temperature has been reached, press the "SET" key under the Heat display or wait a moment and the unit will beep indicating that it has been set. All plate positions will heat to the same settings.
- 5. Press the power button to trun off the unit. **UNPLUG UNIT AFTER EACH USE**.

Potential Hazards













Preventions to reduce exposure to hazards



Vapor Hazard



Eye protection

Protective gloves

No-slip, closed shoes

Read Manual

Heat Hazard

Fire Hazard